


NEW GENERATION CRYOGENIC NITROGEN PLANTS

Your nitrogen, our priority.

The CNP40 is a state-of-the-art liquid nitrogen plant with a production capacity of 40 liters per day. This advanced system also offers a quick startup feature, expediting the process of nitrogen liquefaction while ensuring operational efficiency. Both attributes solidify its standing as a plug-and-liquefy system. Boasting a user-friendly interface and one-button operation, the CNP series liquefiers can be seamlessly integrated into any setup. The necessary electrical connections are all that's required for its fully automated operation, governed by the intelligent PLC controller. The operator's duties are restricted to changing filters and conducting routine checks at maintenance intervals of 8,000 operating hours.



The production of liquid nitrogen is achieved through liquefaction from air, which is then stored in an internal Dewar with capacity of your choice. The availability of liquid nitrogen is ensured at all times, allowing for convenient filling of the dispensing thermos or transfer to external Dewars via a flexible hose with a simple activation. The transfer of liquid nitrogen is independent of the system's operating mode, and the programmable logic controller (PLC) automatically initiates production when the Dewar level drops to 70%. The plant will stop production when the Dewar is full, entering standby mode until liquid nitrogen is transferred.

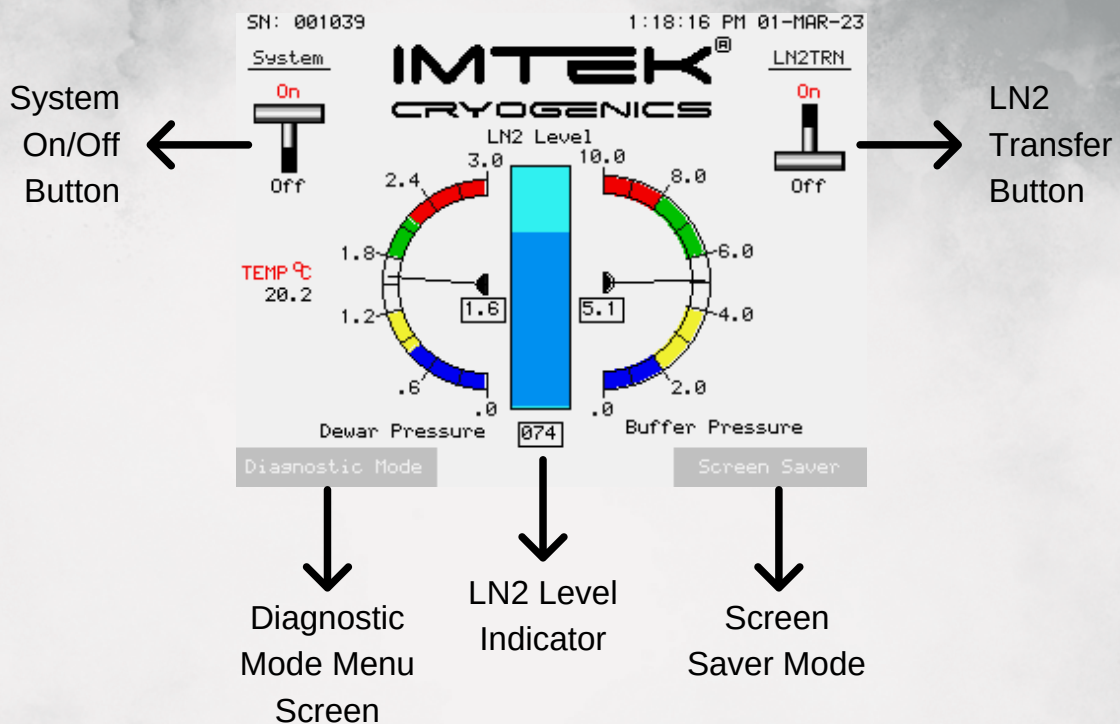

The CNP40 Nitrogen Generation and Storage System's Programmable Logic Controller (PLC) interface includes five essential screens. The Main screen serves as the central control panel, offering start/stop controls, diagnostic mode, and pressure indicators for seamless operation. The Service screen provides direct monitoring and control over key components and features multilingual options. A dedicated Cryo-Service screen ensures customers always have easy access to technical assistance contact information. The Maintenance screen, tailored for system purge and technical interventions, facilitates efficient service operations. Lastly, the Initialization screen aids users during the initial setup and purging, transitioning to the Main screen once the setup is complete. These interfaces seamlessly integrate to provide comprehensive control over the system's operations.



Experience enhanced accessibility and convenience with the CNP40's advanced PLC screen interface. This innovative system not only centralizes control and offers comprehensive monitoring capabilities, but it also enables remote monitoring from your personal computer! Moreover, the Diagnostic's Screen pictured above allows for monitoring of the Air Pressure in the Air Compressor and Buffer Tank as well to ensure safe operation. There is close monitoring of the purity level of the Liquid Nitrogen produced through the O2 Level indicator as well!

The interface persistently exhibits real-time Oxygen and Purity sensor readings, ensuring immediate, accurate understanding of the LN2's quality and enabling quick reactions to any deviations in purity or oxygen levels. Further enhancing its functionality, the PLC interface consistently generates reports encompassing product purity, oxygen levels, system status, and other vital operational parameters. These reports can be configured to be delivered at predetermined intervals or on demand, offering operators a comprehensive insight into the system's performance and the LN2 quality.

02:12:54 PM 24-MAY-23

Buffer AutoPurge		Dewar AutoPurge	
Air Comp OFF	0.0	Air Pressure is LOW	
Buffer Purge OFF	0.0	Dewar Purge OFF	0.0
Circulation Pump OFF		He Comp OFF	
Chiller OFF		CNPTemp is HIGH 0.0	
Restart Level: %	6	Air Comp	00000 h
		He Comp	00000 h


O2 Lvl:
Dew P.:
Level: 0 dec.

English French Spanish

Home CryoService Alarm History Service




Central Line Filter
Water Filter
Particle Filter

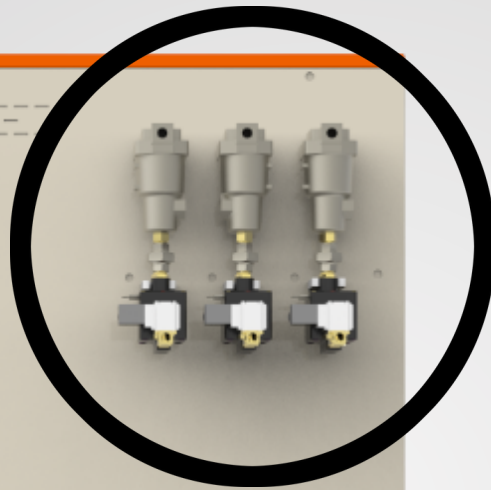


System On/Off
Switch
(Manual)



Left: Water
Inlet/Outlet
Right: Exhaust
Outlets (Water

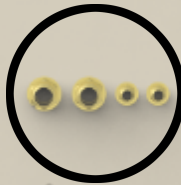

Remote
Control Cable
Connections

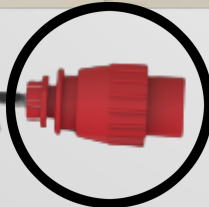

3 Phase
Power Plug



MODEL	CNP40
Production Rate	≥ 40 liter/day (≥48 Lt/day @ 23 °C)
Electrical Options	380 - 415 (±10%) V3 ~ @50 Hz 460 - 480 (±10%) V3 ~ @60 Hz
Power Consumption (Steady State)	3.6 kW @ 50Hz 4.6 kW @ 60Hz
Dimensions (W x L x H)	765 mm x 1340 mm x 1380 mm
Weight	250-300kg
Suggested Installation Area	3m (W) x 3m (L) x 3m (H)
Built-in Air Compressor	Built-In Oil-Free Compressor, ≥ 2 m ³ / hour @7 bar (102 psig)
Cryocooler	GM Type Cryocooler Mounted on Dewar
Compressor	He, 99.995% purity @ 19-19.3bar (275-280 psig), Air Cooled
Req. Cooling Water Flow Rate	4-9 L/min
Req. Chiller Capacity	8 kW
Built-In Nitrogen Generator	
Purity	≥ 99%
Dew Point	up to -40°C
Flow Rate	≥ 1 m ³ / hour
PLC Interface	6" Color Graphic Touch Screen
Dewar Volume	80 Liters (+ Option)
Operating Pressure	1.5 bar
Dewar Level Control	Capacitive Level Sensor
Ambient Temperature Range	+4°C to +40°C
Maximum Altitude	3000 meters
Noise Level	< 65 dB @ 1 meter
Conformities	CE Conformance, ISO 12100:2010, IEC 60204-1, 2006/42/EC, 97/23/EC; ISO9001:2015

CNP is a registered trade mark of Imtek Cryogenics.

We reserve the right to make modifications to our product offerings at any time, without prior notice.